

Алматы (7273)495-231
 Ангартск (3955)60-70-56
 Архангельск (8182)63-90-72
 Астрахань (8512)99-46-04
 Барнаул (3852)73-04-60
 Белгород (4722)40-23-64
 Благовещенск (4162)22-76-07
 Брянск (4832)59-03-52
 Владивосток (423)249-28-31
 Владикавказ (8672)28-90-48
 Владимир (4922)49-43-18
 Волгоград (844)278-03-48
 Вологда (8172)26-41-59
 Воронеж (473)204-51-73
 Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
 Ижевск (3412)26-03-58
 Иркутск (395)279-98-46
 Казань (843)206-01-48
 Калининград (4012)72-03-81
 Калуга (4842)92-23-67
 Кемерово (3842)65-04-62
 Киров (8332)68-02-04
 Коломна (4966)23-41-49
 Кострома (4942)77-07-48
 Краснодар (861)203-40-90
 Красноярск (391)204-63-61
 Курск (4712)77-13-04
 Курган (3522)50-90-47
 Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
 Москва (495)268-04-70
 Мурманск (8152)59-64-93
 Набережные Челны (8552)20-53-41
 Нижний Новгород (831)429-08-12
 Новокузнецк (3843)20-46-81
 Ноябрьск (3496)41-32-12
 Новосибирск (383)227-86-73
 Омск (3812)21-46-40
 Орел (4862)44-53-42
 Оренбург (3532)37-68-04
 Пенза (8412)22-31-16
 Петрозаводск (8142)55-98-37
 Псков (8112)59-10-37
 Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
 Рязань (4912)46-61-64
 Самара (846)206-03-16
 Санкт-Петербург (812)309-46-40
 Саратов (845)249-38-78
 Севастополь (8692)22-31-93
 Саранск (8342)22-96-24
 Симферополь (3652)67-13-56
 Смоленск (4812)29-41-54
 Сочи (862)225-72-31
 Ставрополь (8652)20-65-13
 Сургут (3462)77-98-35
 Сыктывкар (8212)25-95-17
 Тамбов (4752)50-40-97
 Тверь (4822)63-31-35

Тольятти (8482)63-91-07
 Томск (3822)98-41-53
 Тула (4872)33-79-87
 Тюмень (3452)66-21-18
 Ульяновск (8422)24-23-59
 Улан-Удэ (3012)59-97-51
 Уфа (347)229-48-12
 Хабаровск (4212)92-98-04
 Чебоксары (8352)28-53-07
 Челябинск (351)202-03-61
 Череповец (8202)49-02-64
 Чита (3022)38-34-83
 Якутск (4112)23-90-97
 Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://vacuubrand.nt-rt.ru/> || vda@nt-rt.ru

Auxiliary equipment

for vacuum aspiration systems

VacuHandControl VHC^{pro}



Description

The VHC^{pro} is an ergonomic light weight suction hand set, fully autoclavable at 121 °C. It is ready for use with 2.5 m flow through and very flexible silicon tube, adapters for graduated, aspiration and pasteur pipettes, adapter for pipette tips, a wall holder for standby-position (keeps the table top free) and clips with adhesive surface to guide the suction tube into the biosafety cabinet neatly. The mechanism of the VHC^{pro} handset is not in contact with contaminated fluids and enables reliable, leak free working.

Performance features

- light weight and ergonomic for convenient handling
- permanent suction mode for continuous flow
- flow through media tube for leak free working
- autoclavable at 121 °C and 2 bars

Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
VHC ^{pro}	without	-	20688061

VHC^{pro} VHC^{pro}



Description

The stainless steel stand for the VHC^{pro} enables to place the tipped suction handset conveniently and stable during work with cell cultures. It is autoclavable and comes with slip-resistant feet. The stand supports to place the VHC^{pro} also with long, e. g. graduated pipettes, without touching the working surface.

Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
VHC ^{pro}	-	-	20635680

Adapter for pipette tips with BVC ejector



Description

This adapter is designed for pipette tips 2 - 200 µl and has got a tip ejector to remove the used pipette tip

Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
BVC	-	-	20635638

8-channel VHCpro Adapter for Pipette Tips with BVC Ejector



Description

This adapter allows parallel operation of eight pipette tips (size 2 - 200 µl or 5 - 300 µl). It is designed to match the dimensions of standardized 96-well microplates and has got a tip ejector to remove the used pipette tips.

Performance features

- suction power is evenly spread across all pipette tips
- Completely autoclavable
- stand and disinfectant reservoir inclusive

- Low effort for attaching and ejecting the tips

Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
--------------	-----------------------------	------	-----------

BVC

-

-

20635679

Trolley for BVC BVC shuttle

Description

The BVC shuttle is a mobile underframe for all models BVC control and BVC professional. It serves to pull out the BVC easily from under the safety cabinet for cleaning and removal of bottle for change



Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
--------------	-----------------------------	------	-----------

BVC shuttle

-

-

20696880

Adapter for connecting VHC/VHCpro BVC



Description

A set of quick-release connections, with an adapter for connecting a manual manipulator to a container; excellent chemical resistance, equipped with a nozzle to minimize the formation of aerosol and foam, can be used to connect a second manual manipulator. In the disconnected state, it retains tightness and vacuum in the container, autoclavable.

Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
--------------	-----------------------------	------	-----------

BVC

-

-

20635807

Set of connection parts for a VHCpro BVC



Description

Fitting with a nozzle for minimal aerosol and foam formation during use (without a manual manipulator)

Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
BVC	-	-	20699943

Quick-release adapter bottle pump BVC



Description

A set of quick-release connections with an adapter for connecting the container to a BVC or BVC basic system with excellent chemical resistance. In the disconnected state, it retains tightness and vacuum in the container, autoclavable.

Ordering data

Product name	Mains voltage and frequency	Plug	Order no.
BVC	-	-	20635808

BVC PP flask 4L BVC



Description

4L container with high chemical resistance (PP). The front part is designed to reduce the formation of foam and aerosol, a fitting for VHC / VHCpro and a second port for optional connection of an additional VHC / VHCpro adapter, a hydrophobic filter (0.2 microns) to protect the pump and the environment, a hose for connection to a vacuum pump or BVC basic system; fully autoclavable

Ordering data

Product name	Mains voltage and frequency	Plug.	Order no.
--------------	-----------------------------	-------	-----------

BVC glass. capacity 2L BVC**Description**

2L borosilicate glass container is chemically resistant, with an anti-shatter coating, the front part is designed to reduce the formation of foam and aerosol, a fitting for VHCpro and a second port for optional connection of an additional VHCpro adapter holder, a hydrophobic filter (0.2 microns) to protect the pump and the environment, a hose for connection to a vacuum pump or BVC basic system; fully autoclavable.

Ordering data

Product name	Mains voltage and frequency	Plug.	Order no.
BVC	-	-	20635809

Holder for 2L glass bulb BVC**Description**

Holder for 2L glass container for BVC control or BVC professional (included with BVC control G and BVC professional G)

Ordering data

Product name	Mains voltage and frequency	Plug.	Order no.
BVC	-	-	20635839

Spare Hose for VHC/VHCpro VHC/ VHCpro

Silicone hose (minimum order length is 2m)

Ordering data

Product name	Mains voltage and frequency	Plug.	Order no.
--------------	-----------------------------	-------	-----------

VHC / VHC^{pro}

-

-

20636156

Protection filter 0.2 µm 0.2 µm



Description

Hydrophobic 0.2 microns protective filter to protect pump systems; autoclavable up to 20 times

Ordering data

Product name	Mains voltage and frequency	Plug.	Order no.
--------------	-----------------------------	-------	-----------

0.2 µm

-

-

20638266

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Новый Уренгой (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Саранск (8342)22-96-24
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

<https://vacuubrand.nt-rt.ru/> || vda@nt-rt.ru